

# Manual for StillPeaks DHAtOD86 Installation


Thank you for choosing this StillPeaks product. With this manual we'll help you installing.

## Alert: Software and License are separately distributed


Please keep in mind that we distribute the software and the necessary license separately. To get your License first install the software accordingly. Then run the software and follow the License Registration Procedure. After registration we will provide your License-code asap (Please acknowledge 48 hours to process your order/request).

## Download & Install


1. Please download the software from [StillPeaks.com/023020116downloadingpage](http://StillPeaks.com/023020116downloadingpage)  
Be aware to chose the right version for your Windows 10 or Windows 7 computer.
2. Find the Installation-ZIP-file in your downloads and extract it
3. Find the Installation-file in your directory-tree and dubbleclick it

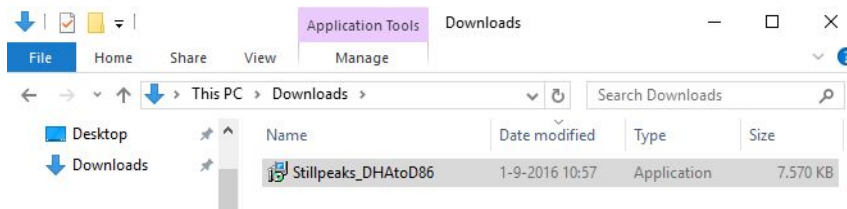


4. Maybe you get a Windows warning about 'Unknown publisher'  
Click '**Run anyway**' or click '**More info**' and then click '**Run anyway**'
5. Start the installation by clicking Next

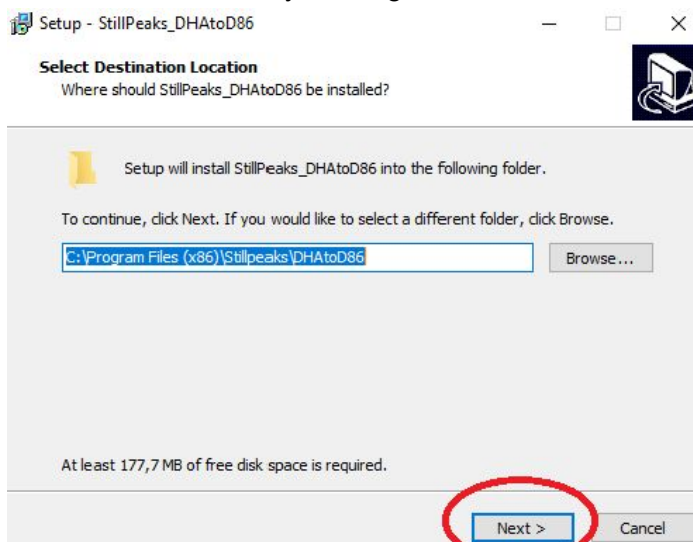


And clicking Next in the following screen:

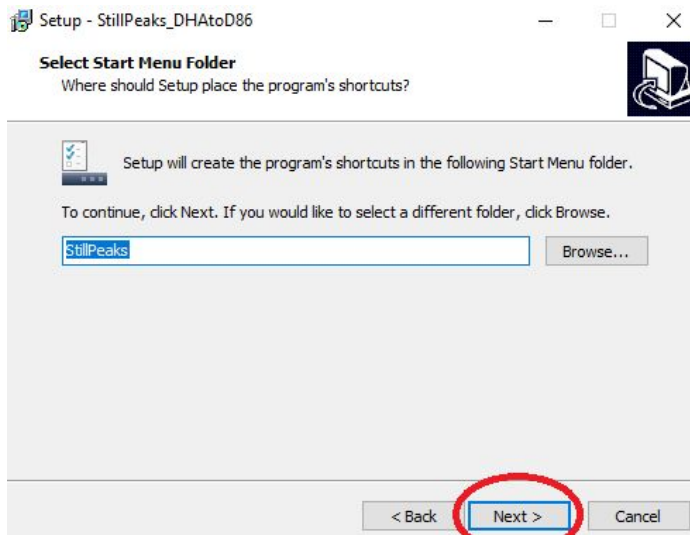

6. Click 'Install'

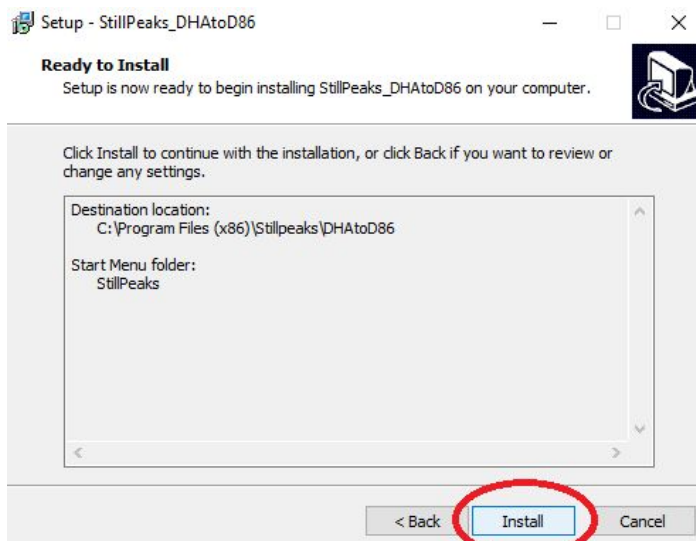


7. Click Finish for completing the installation

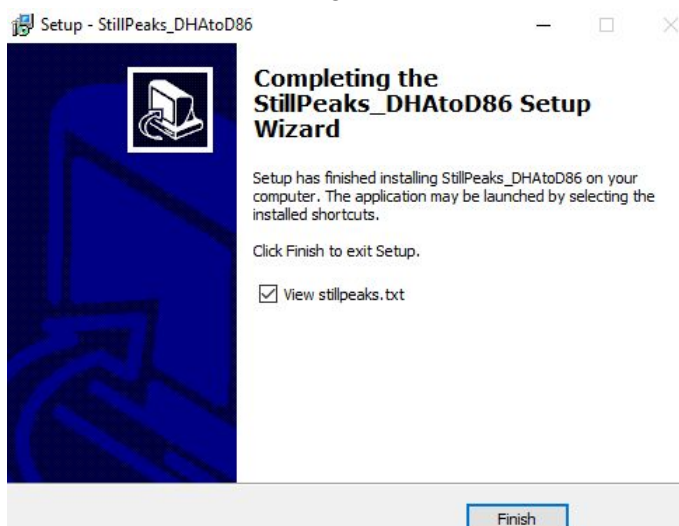



8. Please read the Read-me file called StillPeaks.txt

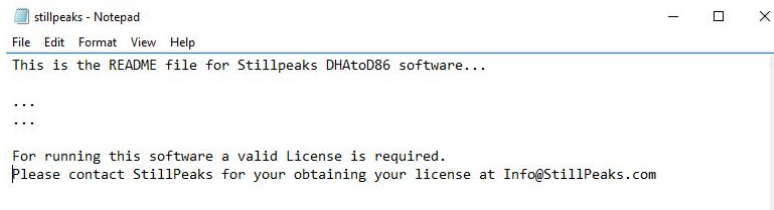

9. Now your StillPeaks Software is installed.

Please check by starting the software via the StillPeaks shortcut

## License

From this point on you require a valid license.

Follow the onscreen instructions and contact us with the appropriate details.

After payment you'll get a valid License-code send. For test purposes we can send you a Free 30 day Trial-License which will provide full access for 30 days.